applied biosystems 7500荧光定量PCR仪

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2.1 输入实验名称 (Experiment Name)

How do you want to identify this experiment?

* Experiment Name:

Barcode (Optional):

User Name (Optional):

Comments (Optional):

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Which instrument are you using to run the experiment?

☒ 7500 (96 Wells) ☐ 7500 Fast (96 Wells)

Set up, run, and analyze an experiment using a 4- or 5-color, 96-well system.

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What type of experiment do you want to set up?

☐ Quantitation - Standard Curve ☐ Quantitation - Relative Standard Curve ☒ Quantitation - Comparative C_T ($\Delta\Delta C_T$)

☐ Melt Curve ☐ Genotyping ☐ Presence/Absence

Use a reference sample and an endogenous control to determine the relative quantity of target nucleic acid sequence in samples.

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Which reagents do you want to use to detect the target sequence?

☒ TaqMan® Reagents ☐ SYBR® Green Reagents ☐ Other

The PCR reactions contain primers designed to amplify the target sequence and a TaqMan® probe designed to detect amplification of the target sequence.

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Which ramp speed do you want to use in the instrument run?

☒ Standard (~ 2 hours to complete a run) ☐ Fast (~ 40 minutes to complete a run)

For optimal results with the standard ramp speed, Applied Biosystems recommends using standard reagents for your PCR reactions.

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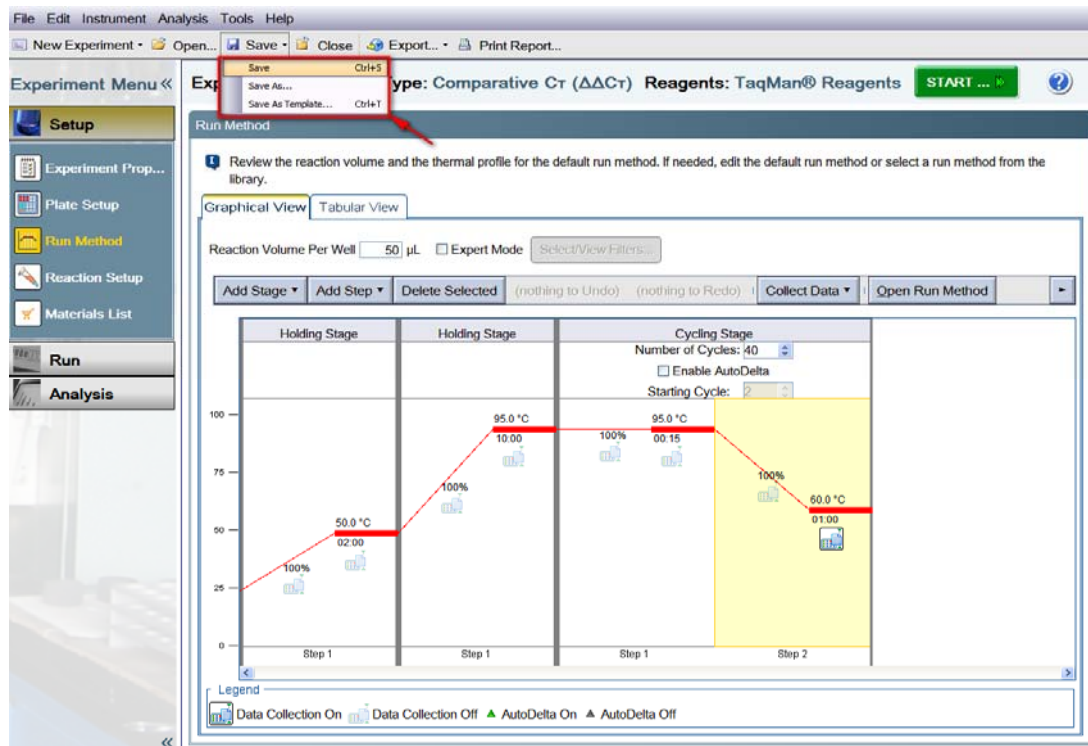
Export Data

Select the type of data to export, select whether to export one file or separate files, then enter export file properties. (Optional) Click "Customize Export" to change the export format and to select fields to export. Click "Start Export" to export your data.

Export Properties | **Customize Export**

1. Select data to export:

- ☐ Sample Setup
- ☒ Results
- ☐ Raw Data
- ☐ Multicomponent Data
- ☐ Amplification Data
- ☐ Melt Region Temperature Data
- ☐ Melt Region Normalized Data
- ☐ Melt Region Derivative Data

2. Select one file or separate files: **One File** Select to export all data in one file or in separate files for each data type.

3. Enter export file properties:

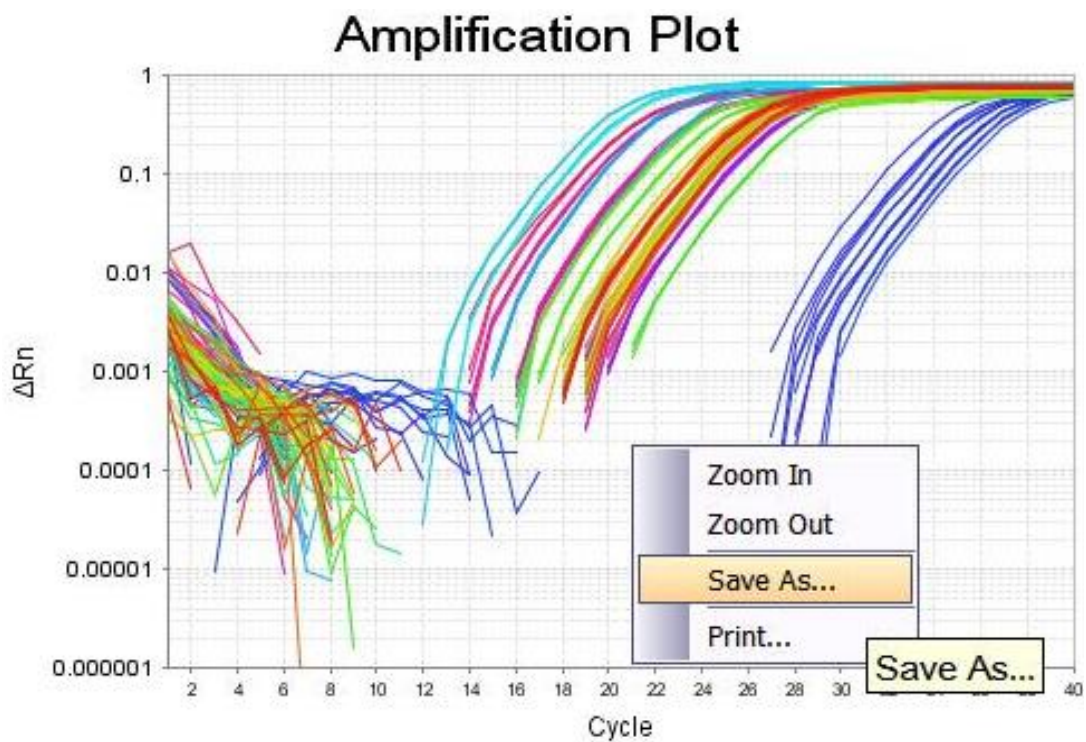
Export File Name: **96-Well Standard Curve Example_data** 更改文件名 File Type: **(* .xls)**

Export File Location: **D:\Applied Biosystems\StepOne Software v2.2\experiments** 更改文件存放位置 **Browse**

☐ Open file(s) when export is complete

☐ Save current settings as the default

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Life Sciences Solutions, Thermo Fisher Scientific

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